



## CELERITY Deluxe SPE Cartridges

Compound: Acetaminophen

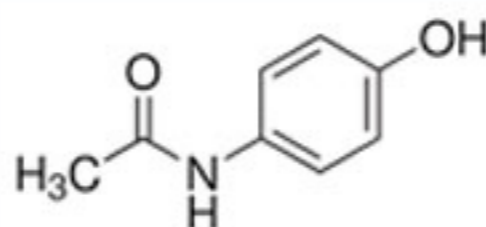


### SPE EXTRACTION METHOD:

(Celerity Deluxe cartridge, 30 mg/1cc)  
(Catalog # SYCY6030-01)

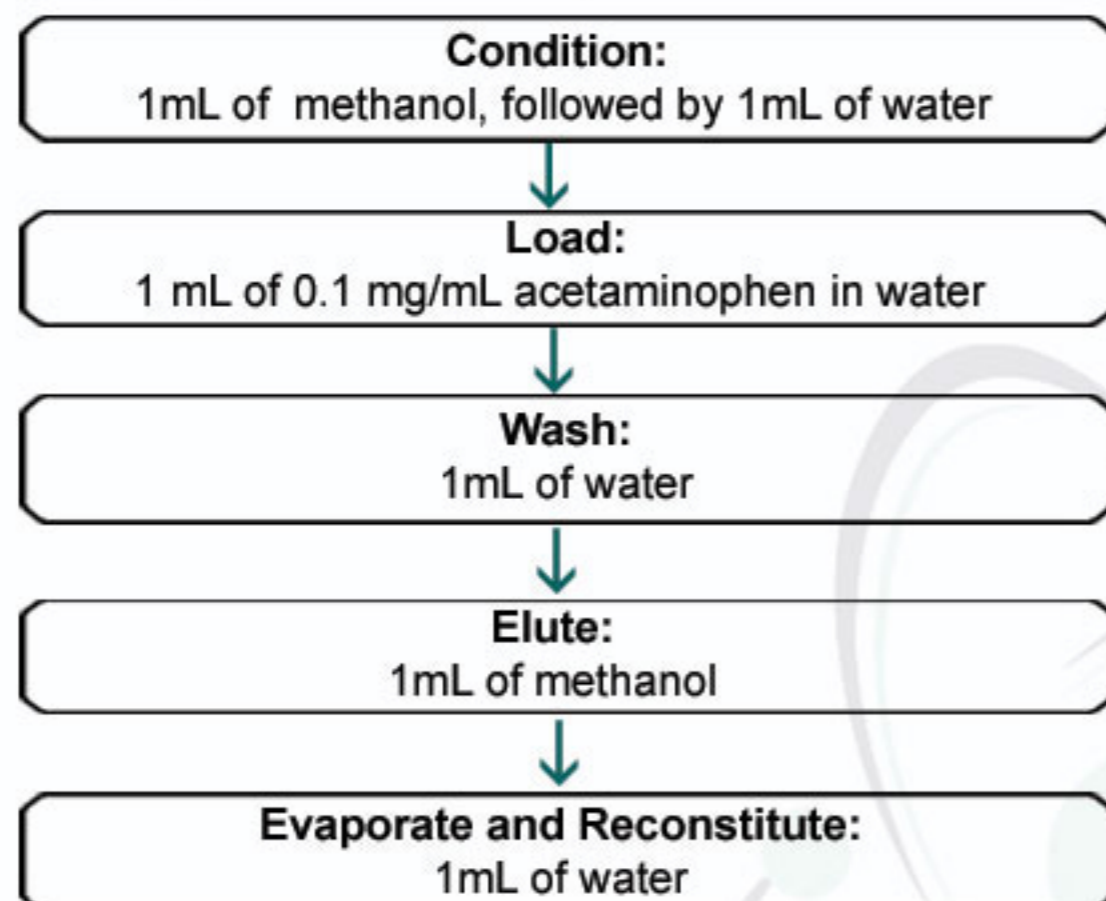

### HPLC METHOD:

**Column:** Orochem Reliasil C18 HPLC Column 4.6x250 mm, 5µm.  
**Part number:** R5HI-121  
**Mobile phase:** Water/acetonitrile/20 mM ammonium acetate 70:20:10  
**Flow rate:** 1mL/min  
**Detection:** UV @220nm  
**Injection volume:** 20µL  
**Retention time:** 4.8 min

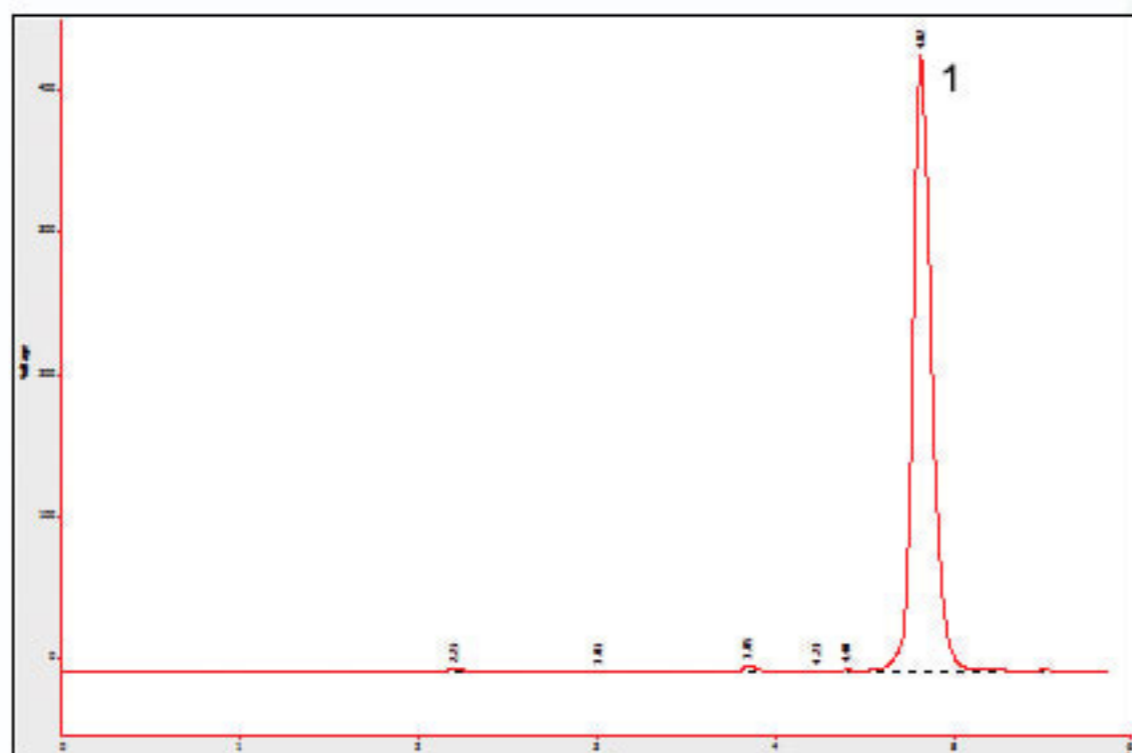

Recovery : 104%