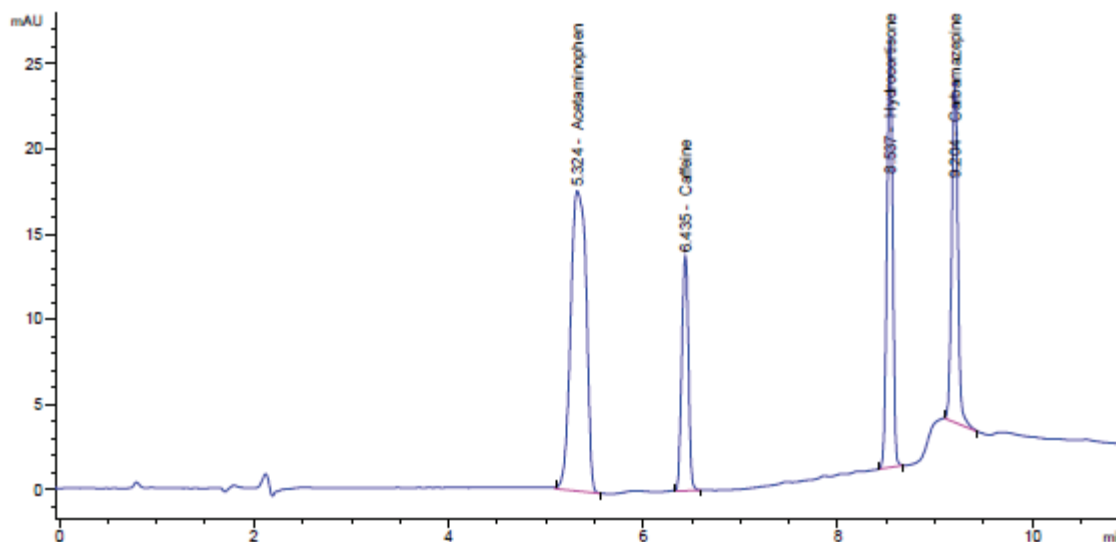
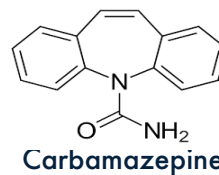
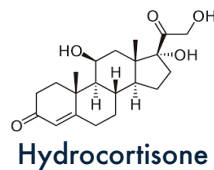
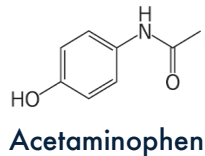
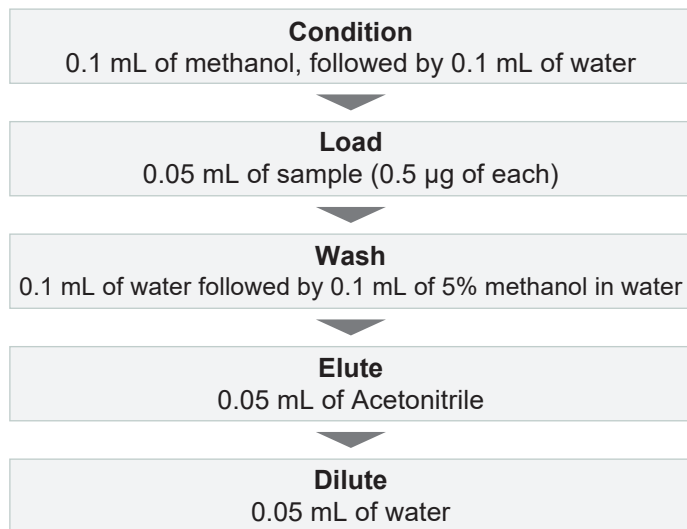


Compounds:

MicroElution Plate Extraction Method:

Panthera Deluxe MicroElute Plate, 3 mg/ well,
 Cat # OC03PA6002


Recovery and precision test results:

	Recovery%	CV(%)
Acetaminophen	106.0	2.6
Caffeine	106.7	2.8
Hydrocortisone	116.2	2.1
Carbamazepine	109.8	2.4

HPLC-UV Method

Column	Orosil C18 HC, 3 µm
Column Size	4.6 mm ID x 150 mm long
Part Number	03FI-023
Temperature	25 °C
Mobile Phase Gradient	A: Water B: ACN Time B (%) 0 10 2 10 8 70 11 70
Injection	10 µL
Flow Rate	1 mL/min
UV Wavelength	254 nm

